Zika Virus NS1 ELISA Quickstart

PREP

- 1. Dilute the 25x wash buffer 1 in 25 with purified lab grade water to make 1x wash buffer
- 2. Reconstitute standards in **BLUE** zero standard as indicated on label and mix well
- 3. Dilute all samples to be tested a minimum of 1 in 10 in **BLUE** zero standard (do not dilute the kit standards)

ASSAY

- 4. Add 100µl of all standards (incl zero standard), samples and any controls to the wells in duplicate
- 5. **Cover plate** with self-adhesive sealing film and incubate with shaking for 2 hours at room temperature
- 6. Aspirate the wells, wash 3 times with 300µl of 1x wash buffer
- 7. Add 100µl PINK Antibody-Biotin reagent to each well, cover plate and incubate with shaking for 1 hour at room temperature
- 8. Aspirate and wash the plate as in step 3
- 9. Add 100µl streptavidin-HRP reagent to each well, cover plate and incubate with shaking for 30 minutes at room temperature
- 10. Aspirate and wash the plate as in step 3, but perform 6 washes
- 11. Add 100µl of TMB substrate to each well, cover plate with a new sealing film and incubate with shaking for 15 minutes at room temperature

.....

12. Add 100µl of stop solution to each well. Read the plate at 450nm, reference at 620nm (605nm-650nm) within 15min of stopping. Use the zero standard as the blank.