

# Prepare UHPLC/HPLC samples faster and minimize leachables with the new glass design

The Whatman™ Mini-UniPrep™ G2 Syringeless filter simplifies UHPLC/HPLC sample preparation over the traditional syringe filter method. The Mini-UniPrep G2 is an all-in-one integrated filter that replaces the syringe, syringe filter, glass vial, cap, and septum (see Fig. 1). After the filtration step Mini-UniPrep G2 can be placed directly into the autosampler in readiness for injecting sample into the UHPLC/HPLC instrument.

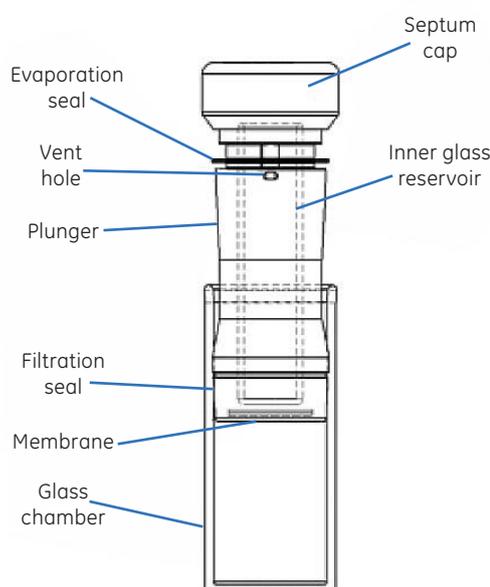


**Fig. 1:** Mini-UniPrep G2 replaces multiple consumables

## Features of the Mini-UniPrep G2:

- Consists of an integral borosilicate glass autosampler vial, plunger with attached filter membrane, and septum/cap
- Filters samples three times faster compared to the traditional syringe filter method
- Glass construction minimizes the risk of leachables contaminating the sample
- Designed to be loaded directly into the autosampler
- Includes visual indication that the sample has been filtered
- Minimizes instrument downtime due to unfiltered samples
- Wide range of membranes with 0.2 and 0.45 µl pore sizes to meet specific sample filtration requirements

The Mini-UniPrep G2 includes an integral borosilicate glass vial housed within the plunger (see Fig. 2) and a borosilicate glass chamber for holding the unfiltered liquid. During the filtration step, the plunger is compressed into the glass chamber containing the unfiltered liquid. As the plunger travels downward, liquid flows through the filtration membrane to the top of the plunger and drops into the glass collection vial housed within the plunger (see Fig. 3). Therefore, the sample only contacts plastic for a very short period of time i.e. only while the plunger is being compressed through the unfiltered liquid. Once compressed, the Mini-UniPrep G2 is ready to be loaded directly on to the autosampler.

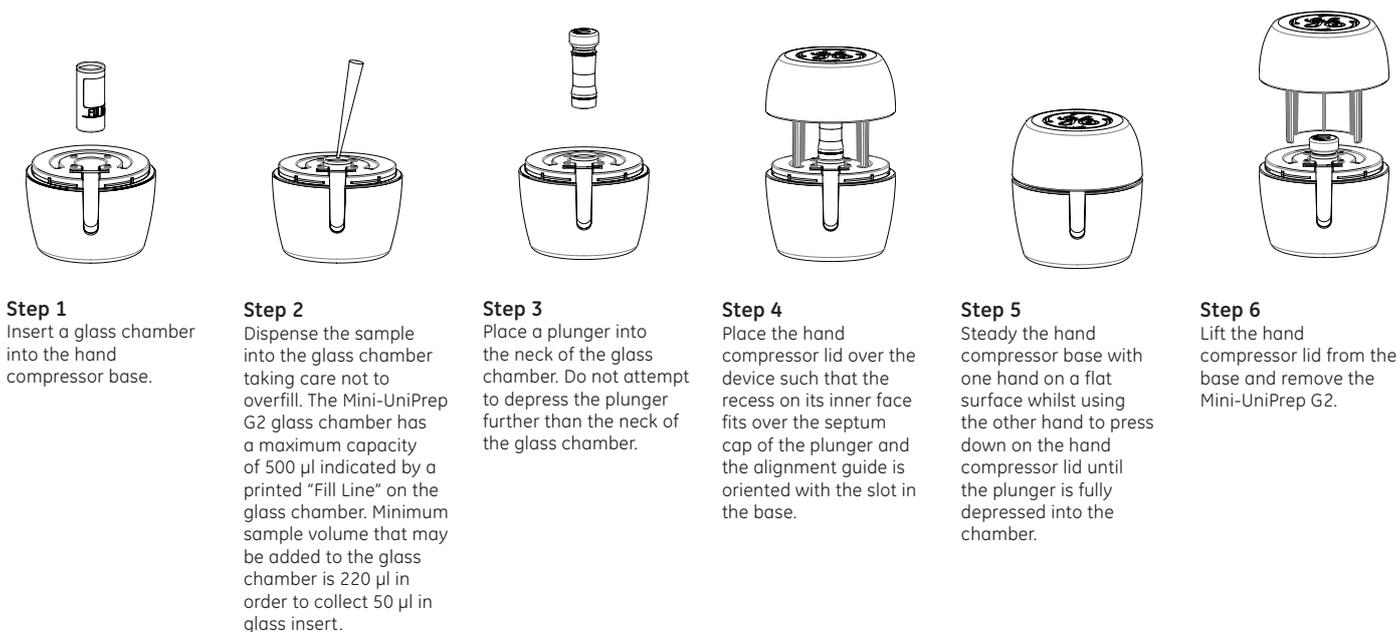


**Fig. 2:** Mini-UniPrep G2

Mini-UniPrep G2 comes in a number of configurations including slit septa for those autosamplers and robotic systems that require a slit septa. It is also available in amber for light sensitive samples. Amber Mini-UniPrep meets the requirements of the US Pharmacopeia (USP) and European Pharmacopeia (EP) for light transmission.



**Fig. 3:** How to use the Mini-UniPrep G2



## Technical Specifications

<b>Dimensions</b>	Once compressed, equivalent in size to 12 mm x 32 mm vial
<b>Materials of Construction</b>	Chamber: Borosilicate Glass Plunger Outer Housing: Polypropylene Plunger Inner Storage Vial: Borosilicate Glass Filter Media: As specified Septa: Silicone with PTFE Liner Cap: Polypropylene
<b>Filtering Capacity:</b>	Chamber (Unfiltered Sample): 500 µl Inner Storage Vial (Filtered Sample) : 330 µl Recommended Minimum Filtering Volume: 220 µl placed in the Chamber to obtain 50 µl in Inner Storage Vial
<b>Nominal Force Needed to Compress</b>	Approx. 25 lbs (11.3 Kg)
<b>Autosampler Compatibility</b>	Any autosampler that accommodates standard 12 mm x 32 mm profile vials (Needle height of the autosampler may need adjusting)

For local office contact information, please visit [www.gelifesciences.com/contact](http://www.gelifesciences.com/contact)

GE Healthcare UK Limited  
Amersham Place  
Little Chalfont  
Buckinghamshire, HP7 9NA  
UK



GE, imagination at work and GE monogram are trademarks of General Electric Company.

Mini UniPrep and Whatman are trademarks of GE Healthcare companies.

©2012 General Electric Company—All rights reserved.

First published August 2012.

All goods and services are sold subject to the terms and conditions of sale of the company within GE Healthcare which supplies them. A copy of these terms and conditions is available on request. Contact your local GE Healthcare representative for the most current information.

GE Healthcare Bio-Sciences Corp.  
800 Centennial Avenue, P.O. Box 1327  
Piscataway, NJ 08855-1327  
USA

GE Healthcare Bio-Sciences AB  
Björkgatan 30  
751 84 Uppsala  
Sweden

GE Healthcare Europe, GmbH  
Munzinger Strasse 5  
D-79111 Freiburg  
Germany

GE Healthcare Japan Corporation  
Sanken Bldg., 3-25-1, Hyakunincho  
Shinjuku-ku, Tokyo 169-0073  
Japan