

219 North Arkansas Street Rogers, AR 72756

Phone: (479) 636-4361; 800-643-3426 Fax: (479) 636-3562

Email: biosales@pel-freez.com
Website: www.pelfreez-bio.com

PRODUCT DATA SHEET

3-4 WEEK RABBIT COMPLEMENT

Product Codes	Pack size
31061-0	1 mL
31061-1	10 mL
31061-3	100 mL

Description:

Complement is produced under controlled conditions from large pools of New Zealand White rabbits that are 21-28 days old. This provides a product that has low heterophile antibody activity for many applications, including use in bactericidal and opsonization assays.

Physical State:

Frozen liquid

Testing:

Hemolytic: Titer ≥ 1:16 Endotoxin: ≤ 1.0 EU/mL

Sterility: Passed, 14 day sterility testing

Mycoplasma: Negative

Viral: Negative

Packaging, shipping/storage:

Packaging

1 mL: Glass vial; 10 mL and 100 mL: PETG bottle

Storage Temperature

-70 °C or below

Shipping Conditions

Dry ice

Expiration

Product quality is guaranteed to meet Pel-Freez Biologicals' specifications for 1 year from the date of receipt by the customer as long as the product is stored in accordance with the indicated storage conditions.

Application Notes:

Suitable for functional antibody assays (SBA-Serum Bactericidal Assay, MOPA/OPA- Multiplexed /Opsonophagocytic Killing Assay) for vaccine evaluation. Complement components are extremely heat-sensitive; for best results, thaw complement on ice or in a cold water bath (ideally thaw bottled complement under a stream of cold water 20°C); never thaw using warm or hot water. Single use aliquots may be made at this time and frozen at -70°C or below. Avoid multiple freeze-thaw cycles. When a working aliquot is thawed, do not re-freeze remaining complement, but discard it.

P/N: 74012 Rev 02 Page 1 of 2



219 North Arkansas Street Rogers, AR 72756

Phone: (479) 636-4361; 800-643-3426 Fax: (479) 636-3562

Email: biosales@pel-freez.com Website: www.pelfreez-bio.com

References:

Hirve S, Bavdekar A, Pandit A, Juvekar S, Patil M, Preziosi MP, Tang Y, Marchetti E, Martellet L, Findlow H, Elie C, Parulekar V, Plikaytis B, Borrow R, Carlone G, Kulkarni PS, Goel A, Suresh K, Beri S, Kapre S, Jadhav S, Preaud JM, Viviani S, LaForce FM. Immunogenicity and safety of a new meningococcal A conjugate vaccine in Indian children aged 2-10 years: a phase II/III double-blind randomized controlled trial. Vaccine. 2012 Oct 5;30(45):6456-60.

Romero-Steiner, S., C. E. Frasch, G. Carlone, R. A. Fleck, D. Goldblatt, and M. H. Nahm. 2006. Use of opsonophagocytosis for serological evaluation of pneumococcal vaccines. Clin. Vaccine Immunol. 13:165-169.

Romero-Steiner, S., D. LiButti, L.B. Pais, J. Dykes, P. Anderson, J.C. Whitin, H.L. Keyserling, and G.M. Carlone. 1997. Standardization of an opsonophagocytic assay for the measurement of functional antibody activity against *Streptococcus pneumoniae* using differentiated HL-60 cells. Clin. Diagn. Lab. Immunol. 4:415-422.

Burton R.L., Nahm M.H.. Clin Vaccine Immunol. 2012 Jun;19(6):835-41. Development of a fourfold multiplexed opsonophagocytosis assay for pneumococcal antibodies against additional serotypes and discovery of serological subtypes in Streptococcus pneumoniae serotype 20.

P/N: 74012 Rev 02 Page 2 of 2